## LAUTSTÄRKEREGELUNG

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## TIMER FUNCTIONS & DESCRIPTION

### M1 button (1)

Press the button to set the first storage time. Press and hold the button for 2 seconds to access the countdown time setting.

### M2 button (2)

Press the button to set the second memory time. Press and hold the button for 2 seconds to access the countdown time setting.

### M3 button (3)

Press the button to set the second storage time. Press and hold the button for 2 seconds to access the countdown time setting.

### MIN button (4)

Press the button to change the minutes in setting mode.

### SEC button (5)

Press the button to change the seconds in setting mode.

### S/S button (6)

Press the S/S button to start and stop the timer function.

### Mute / Lo/Hi button (7)

Slide button to change the volume.

Wall bracket (8), magnet (9), stand (10), battery cover (11)

## LCD-DISPLAY

Time window (1), minute (2), second (3), count-up indicator (4), count-down indicator (5)

## VOLUME CONTROL

You can control the volume of the signal using the side slide button (7). You can choose between loud (Hi), quiet (Lo) and mute (Mute).

## MOUNTING OPTIONS FOR THE TIMER

Use the wall bracket (8), the magnet (9) or the stand (10) on the back of the appliance for positioning. Mounting material for wall mounting is not included in the scope of delivery.

## COUNTDOWN MEMORY SETTING

1. Press and hold the M1 button (M2 or M3) for 2 seconds to access the timer time setting.
2. Press the MIN button to set the minutes.
3. Press the SEC button to set the seconds.
4. Press the M1 button (M2 or M3) again to confirm and save the setting. This also happens automatically after 30 seconds in which no button is pressed.
5. Repeat steps 1-4 to change the saved data.
6. Press and hold the MIN and SEC buttons simultaneously to set the count data to 0.

Count time mode

Count down memory setting

## USING THE COUNTDOWN FUNCTION

1. Set the desired countdown time.
2. Press the M1 button (M2 or M3) to select the saved time. Now press the S/S button to start or stop the countdown function. While the countdown is running, the light flashes green.
3. The alarm sounds as soon as the countdown time reaches 0. When the time has elapsed, the light flashes red. Press the S/S button to acknowledge the alarm and repeat the saved time. The alarm tone stops automatically after one minute.
4. The timer automatically switches to the count-up function after the count-down time has elapsed.

## USING THE COUNT-UP FUNCTION

If 00<sub>M</sub>00<sub>S</sub> appears on the display, press the S/S button to start the count-up function. Press the S/S button again to stop the timer. This process can be repeated as often as required. The maximum display is 99 minutes 59 seconds.

## DECLARATION OF CONFORMITY

A declaration of conformity and the digital version of these operating instructions can be found here: <http://www.thgeyer.com/declaration-of-conformity-timer>

## BATTERY RETURN

Batteries must not be disposed of with household waste. You are legally obliged to return batteries to suitable collection centres. Batteries containing hazardous substances are labelled with the 'crossed-out dustbin' symbol and one of the chemical symbols Cd (cadmium), Hg (mercury) or Pb (lead).

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