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T1/T2-Taste (1/2)

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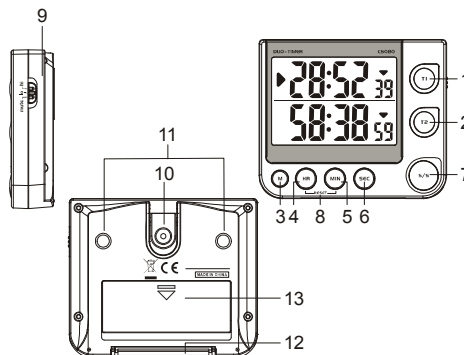
TIMER FUNCTIONS & DESCRIPTION

T1/T2 button (1/2)

Press one of the buttons to set the respective storage time.

M button (3)

Press and hold the button for 2 seconds in the respective memory time to access the countdown time setting. Press again to save the entered value.

HR button (4)

Press the button to change the hours in setting mode.

MIN button (5)

Press the button to change the minutes in setting mode.

SEC button (6)

Press the button to change the seconds in setting mode.

S/S button (7)

Press the S/S button to start and stop the timer function.

Reset (8)

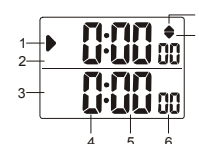
Press the MIN and HR buttons simultaneously to delete all values. In the event of a malfunction, the timer must be reset completely.

Mute / Lo/Hi button (9)

Slide button to change the volume.

Wall bracket (10), magnet (11), stand (12), battery cover (13)

LCD-DISPLAY

Indicator for activated timer time (1), display for timer 1 (2), display for timer 2 (3), hours (4), minutes (5), seconds (6), count-down indicator (7), count-up indicator (8)

VOLUME CONTROL

You can control the volume of the signal using the side slide button (9). You can choose between loud (Hi), quiet (Lo) and mute (Mute).

MOUNTING OPTIONS FOR THE TIMER

Use the wall bracket (10), the magnet (11) or the stand (12) on the back of the appliance for positioning. Mounting material for wall mounting is not included in the scope of delivery.

Count time mode

Count down memory setting

COUNTDOWN MEMORY SETTING

1. In the Timer 1 or Timer 2 display, press and hold the M button for 2 seconds until the numbers flash to access the memory setting.
2. Press the HR, MIN or SEC button to set the hours, minutes or seconds.
3. Press the M button again to confirm and save the setting. This also happens automatically after 30 seconds in which no button is pressed.
4. Repeat steps 1-3 to change the saved data.
5. Press and hold the HR and MIN buttons simultaneously (reset) to reset the count data to 0.

USING THE COUNTDOWN FUNCTION

1. Set the desired countdown time.
2. Press the T1 or T2 button to select the saved time. Now press the S/S button to start or stop the countdown function. While the countdown is running, the light flashes green.
3. The alarm sounds as soon as the countdown time reaches 0. When the time has elapsed, the light flashes red. Press the S/S button to acknowledge the alarm and repeat the saved time. The alarm tone stops automatically after one minute.
4. The timer automatically switches to the count-up function after the count-down time has elapsed.

USING THE COUNT-UP FUNCTION

If 0:00₀₀ appears on the display, press the S/S button to start the count-up function. Press the S/S button again to stop the function. This process can be repeated as often as required. The maximum display is 99 hours 59 minutes and 59 seconds.

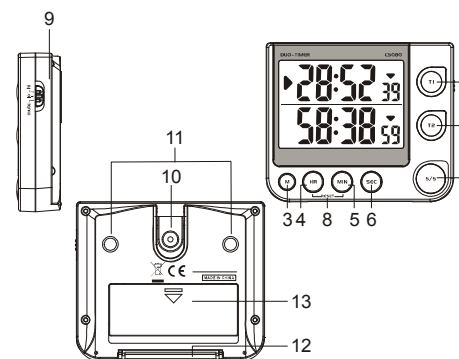
DECLARATION OF CONFORMITY

A declaration of conformity and the digital version of these operating instructions can be found here: <http://www.thgeyer.com/declaration-of-conformity-timer>

BATTERY RETURN

Batteries must not be disposed of with household waste. You are legally obliged to return batteries to suitable collection centres. Batteries containing hazardous substances are labelled with the 'crossed-out dustbin' symbol and one of the chemical symbols Cd (cadmium), Hg (mercury) or Pb (lead).

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